PH ELECTRODES



•	-		_					_			_
GE 100	GE 101	GE 104	GE 108	GE 114	GE 117	GE 120	GE 125	GE 126	GE 151	GE 171	GE 173
0 14 pH 0 80 °C	2 11 pH 0 60 °C	0 14 pH 0 80 °C	0 14 pH 0 80 °C	0 14 pH 0 60 °C	0 14 pH 0 80 °C	0 14 pH 0 60 °C	0 14 pH 0 70 °C	0 14 pH -5 +80 °C	0 14 pH 0 80 °C	0 14 pH 0 140 °C	0 14 pH 0 80 °C
>100 µS/cm	>100 µS/cm	>20 µS/cm	$>100\mu\text{S/cm}$	>200 µS/cm	>100 µS/cm	>200 µS/cm	>200 µS/cm	$>100~\mu S/cm$	$>100\mu\text{S/cm}$	>100 µS/cm	>50 μS/cm
no	no	no	no	no	integr. Pt1000 4 mm banana	no	integr. Pt1000 4 mm banana	no	no	no	no
no	no	no	no	optional	no	no	ja	no	no	no	no
no	no	no	6 bar	no	6 bar	no	1 bar	5.5 bar	no	10 bar	6 bar
1 m 1)	1 m 1)	1 m 1)	2 m 1)	1 m	2 m ²⁾	1 m	2 m	5 m	1 m 1)	ohne	1 m 1)
3 mol/l KCl	3 mol/l KCl	3 mol/l KCl	gel electrolyte	gel electrolyte	gel electrolyte	gel electrolyte	gel electrolyte	gel electrolyte	3 mol/l KCl	gel electrolyte	gel electrolyte
2 x ceramic	2 x ceramic	moving joint	2 x ceramic	1 x Pellon	2 x ceramic	2 x ceramic	1 x ceramic	2 x ceramic	1 x ceramic	2 x ceramic	joint
without	without	without	PG 13.5	without	PG 13.5	without	without	1/2" NPT	without	PG 13.5	PG 13.5
tyril, Ø 12 mm x 120 mm	glass, Ø 12 or 6 mm x 120 mm	glass, Ø 12 mm x 120 mm	PSU, Ø 12 mm x 120 mm	epoxide, Ø 12 mm x 120 mm	PSU, Ø 12 mm x 120 mm	PVC, Ø 22 mm x 110 mm	epoxide, Ø 12 mm x 120 mm	ABS Ø 26.4 mm x 147 mm	glass, Ø 12 mm x 120 mm	glass, Ø 12 mm x 120 mm	glass, Ø 12 mm x 120 mm
universal electrode	tip Ø 6 mm, small sample volume	for low-ion media	low- maintenance	Low-cost low- maintenance	temperature compen- sated	insertion electrode, blade Ø 13 mm x 60 mm	submersible, water-proof IP67 (also BNC-plug)	extremely low- maintenance	chemicals- resistant glass shaft	for extreme conditions, sterilizable, autoclavable	for process chemistry, bio-che- mistry, alkali- resistant
600704	600693	602063	600713	604701	600730	600698	600731	610987	600727	-	600735
-	-	-	606089	-	-	-	-		-	606375	606572
	0 14 pH 0 80 °C >100 μS/cm no no 1 m 1) 3 mol/l KCl 2 x ceramic without tyril, Ø 12 mm x 120 mm universal electrode	0 14 pH 2 11 pH 0 80 °C 0 60 °C >100 μS/cm >100 μS/cm no no no no 1 m 1) 1 m 1) 3 mol/l KCl 3 mol/l KCl 2 x ceramic without tyril, Ø 12 mm x 120 mm universal electrode volume 2 11 pH 0 12 m ly H 100 μS/cm 1 m 10	0 14 pH 0 14 pH 0 14 pH 0 80 °C 0 60 °C 0 80 °C >100 μS/cm >100 μS/cm >20 μS/cm no no no no no no	0 14 pH 2 11 pH 0 14 pH 10 10 pK 0 log	0 14 pH 2 11 pH 0 14 pH 0 16 v C 10 color 0 10 color 0 10 color 0 color 0 100 μS/cm >200 μS/c	0 14 pH 2 11 pH 0 14 pH 0 16 °C 0 80 °C	0 14 pH 2 11 pH 0 14 pH 10 pH 1000	014 pH 016 °C 080 °C 0	014 pH 0	014 pH 080 °C 060 °C 014 pH 014 pH 014 pH 014 pH 080 °C 080 °C 080 °C 080 °C 080 °C 080 °C 060 °C 070 °C 070 °C 014 pH 014 pH 080 °C 070 °C 070 °C 080 °C 0	014 pH 080 °C 060 °C 080 °C 080 °C 080 °C 060 °C 080 °C 0

^{*)} Note: cable GEAK-2S7-BNC or GEAK-5S7-BNC is needed for connection S7, for devices with cinch connection adapter GAD 1 BNC is necessary. Electrodes are consumption objects. Lifetime under careful treatment: >2 years; warranty: 12 months

Options:

Longer cable for 1) 2)

(available cable lengths: up to 5 m)

Special designs

(electrodes with thread, other lengths, special applications etc.)

Accessories and spare parts:

Kabel-BNCM/BNCF

Art. no. 606158

Extension cables for electrodes with BNC connector, Cable length: 3 m



S7 connection at shaft

Diaphragm:

The diaphragm makes the electric connection between reference system and sample. Additionally it should prevent the spoiling of the reference electrolyte by the measured medium.

Ceramic diaphragm

Porous ceramic rods ensure low leak rates

Application:

General applications in clean till lightly soiled media.

Joint / movable joint

The roughened surface between the cut glass of the electrode and a cut glass sleeve permits a electrolyte flow of several ml/h.

Application:

low-ion or heavily soiled samples

Pellon diaphragm

A permeable diaphragm made of Pellon texture is used for fast response times and stable measuring values

Application:

Clean till lightly soiled media.

Reference electrolyte:

The reference electrolyte offers a constant voltage of the reference system and makes the electrical connection between sample and reference electrode.

Liquid electrolyte

Mainly 3 mol/I KCl is used. Liquid electrolytes offer fast response times in general and can be replaced if contaminated.

Gel electrolyte

ceramic rod

glass sleeve

The electrolyte is solidified for low-maintenance electrodes able to measure irrespective to its position. Under normal measurement conditions no noticeable electrolyte leakage is observable.

Electrodes with S7 connection:

The electrodes are offered with an S7 industrial screw plug fitted, also known as industrial-S8 Plug head. In contrast to S7 lab plug head this one is for direct installation in fittings with PG 13.5 suitable thread.



APPLICATION AREAS: ELECTRODES

APPLICATION	GE 100	GE 101	GE 104	GE 108	GE 114	GE 117	GE 120	GE 125	GE 151	GE 171	GE 173	GE 126	GR 105	GR 175
Sewage													Ĭ	
Aquarium water	•		•	•	•	•			•			•	•	•
Soil testing		•												
Emulsions		•	•											
On-site measurements				•	•	•		•					•	
Fish farming	•		•	•	•	•		•	•			•	•	•
Galvanic baths											•			•
Beverages								•	•		•		•	•
Low-ion media			•								•			
Cosmetics			•											
Food sample		•					•							
Sea water	•	•	•	•	•	•	•	•	•	•	•	•	•	•
Online measuring										•	•			•
Process chemistry									•	•	•			•
Swimming pool water	•			•	•	•		•			•	•	•	•
Suspensions		•	•											•
Drinking water	•		•	•	•	•		•			•	•	•	•
Water-based lacquers			•								•			

Note: The set information are to provide general recommendations. It needs to be checked, which electrodes for each area of application are suitable.

ORP ELECTRODES



Art. no. 607798

ORP / redox electrode with BNC connection



GR 175-BNC

Art. no. 607801

ORP electrode with BNC connection

GR 175-S7

Art. no. 607802

ORP electrode incl. S7 connector-without connecting cable *

Electrodes are consumption objects. Lifetime under careful treatment: >2 years; warranty: 12 months

Specifications:	GR 105	GR 175					
Measuring unit:	ORP						
Measuring range:	±2000 mV, 0 80 °C						
Conductivity:	>100 µS/cm						
Temperature measurement:	no						
Water-proof:	no						
Pressure resistant:	no	6 bar					
Cable:	1 m 1)	without/1 m					
Electrolyte:	3 mol/I KCL	gel electrolyte					
Diaphragm:	2 x ceramic	1 x ceramic					
Metal electrode:	Platin dome Ø 5 mm						
Thread:	without	PG 13.5					
Electrode shaft:	tvril. Ø 12 mm	glass. Ø 12 mm					

x 120 mm x 120 mm Minimal depth of immersion: 15 mm

Options:

Longer cable for 1) 2)

Scope of supply:

(available cable lengths: up to 5 m)

Accessories:

GRP 100

Art. no. 601424

Redox test solution 220 mV, 100 ml



ORP electrode, manual

ELECTRODES - ACCESSORIES

GPH 4,0 / 5 pH buffer capsules (5 pieces), pH 4.0	Art. no. 60261
GPH 4,0 / 10 pH buffer capsules (10 pieces), pH 4.0	Art. no. 60261
GPH 7,0 / 5 pH buffer capsules (5 pieces), pH 7.0	Art. no. 60261
GPH 7,0 / 10 pH buffer capsules (10 pieces), pH 7.0	Art. no. 60261
GPH 10,0 / 5 pH buffer capsules (5 pieces), pH 10.0	Art. no. 60261
GPH 10,0 / 10 pH buffer capsules (10 pieces), pH 10.0	Art. no. 60261
GPH 12,0 / 5 pH buffer capsules (5 pieces), pH 12.0	Art. no. 60262
GPH 12,0 / 10 pH buffer capsules (10 pieces), pH 12.0	Art. no. 60262
All buffer capsules are traceable to NIST have ± 0.02 pH at 25 °C.	standards and
PHL 4 pH buffer solution, ready to use (pH 4.0	Art. no. 60136 1 / 25 °C), 250 ml
PHL 7 pH buffer solution, ready to use (pH 7.0)	Art. no. 60137 0 / 25 °C), 250 ml
PHL 10 pH buffer solution, ready to use (pH 10.0	Art. no. 60137 1 / 25 °C), 250 ml
PHL 4-1000 pH buffer solution, ready to use, (pH 4.01	Art. no. 41502 / 25 °C), 1000 ml
PHL 7-1000 pH buffer solution, ready to use, (pH 7.00	Art. no. 41503 / 25 °C), 1000 ml
PHL 10-1000 pH buffer solution, ready to use, (pH 10.0)	Art. no. 41503 1 / 25 °C), 1000 ml
KCL 3 M Art. no. 602477 3 for refilling and storage (fill into protective) with 3 mol KCl electrolyte, injection bot	1.7
CaCl 1000 ml, solution for measuring the pH	Art. no. 60325 value of soil
GRL 100 HCL/Pepsin cleaning solution, 100 ml	Art. no. 60142

GEAK-5S7-BNC

Art. no. 601998

Adapter cable pH S7-BNC, 5 m

VD120

Art. no. 601380

injection aid for injection electrode GE101

GAD 1 BNC

Art. no. 601382

Adapter for the plug-in of electrodes with Cinch plug to devices with BNC socket.

GPF 100

Art. no. 601417

Plastic bottle with wide neck, 100 ml

GAK 1400

Art. no. 603523

pH Working and calibration set: GPH 4,0, GPH 7,0, GPH 10,0 (5 capsules of each type); 3 x GPF 100; 1 x KCL 3M; 1 x GRL 100

GWA1Z

Art. no. 602914

pressure accessories, plastic adapters PG13.5 to G1"

PG 13.5

Art. no. 603205

Thread Adapter, pluggable, Pg 13,5 for sensors with shaft Ø 12 mm



Art. no. 606171

pressure accessories, plastic adapters from PG11 external thread to PG 13.5 internal thread incl. sealing and PG11 counter nut, material: polyamide, fiber glass reinforced, O-ring: NBR, temperature range: -10 ... +100 °C

HD-22-3

Art. no. 700040 freely positionable laboratory sensor holding arm for sensors Ø12mm

SDW 500 Deionized water in refillable 500 ml bottle

^{*)} Note: cable GEAK-2S7-BNC or GEAK-5S7-BNC is needed for connection S7, for devices with cinch connection adapter GAD 1 BNC is necessary.