



HD 2156.1, HD 2156.2
pH METER - CONDUCTIVITY METER - THERMOMETER

The **HD2156.1** and **HD2156.2** are portable instruments with a large LCD display. They measure pH, mV, redox potential (ORP), conductivity, liquid resistivity, total dissolved solids (TDS) and salinity using combined 4-ring and 2-ring conductivity/temperature probes. Temperature only is measured by Pt100 or Pt1000 immersion, penetration, contact or air probes.

The pH electrode calibration, as well as manual, can be carried out on one, two or three points and the calibration sequence can be chosen from a list of 13 buffers.

The calibration of the conductivity probe can be performed automatically in one or more of the 147µS/cm, 1413 µS/cm, 12880 µS/cm or 111800 µS/cm solutions.

The HD2156.2 instrument is a **datalogger**. It stores up to 20,000 sets of three measurements composed of pH or mV, conductivity or resistivity or TDS or salinity and temperature: these data can be transferred to a PC from the instrument connected via the RS232C or USB 2.0 serial ports. The storing interval, printing, and baud rate can be configured by the menu.

Both models are fitted with an RS232C serial port and can transfer the acquired measurements to a PC or to a portable printer in real time.

The *Max*, *Min* and *Avg* function calculates the maximum, minimum or average values.

Other functions include: the Auto-HOLD function and the automatic turning off which can also be excluded.

The instruments have IP66 protection degree.

INSTRUMENT TECHNICAL CHARACTERISTICS

Measured quantities: pH, mV, χ , Ω , TDS, NaCl, °C, °F

Instrument

| | |
|---|---|
| Dimensions (Length x Width x Height) | 185x90x40mm |
| Weight | 470g (complete with batteries) |
| Materials | ABS, rubber |
| Display | 2x4½ digits plus symbols Visible area: 52x42mm |

Operating conditions

| | |
|---------------------------|--------------------------------|
| Working temperature | -5...50°C |
| Storage temperature | -25...65°C |
| Working relative humidity | 0...90%RH without condensation |
| Protection degree | IP66 |

Power

| | |
|------------------------------------|---|
| Batteries | 4 1.5V type AA batteries |
| Autonomy | 200 hours with 1800mAh alkaline batteries |
| Power absorbed with instrument off | 20µA |
| Mains-supply unit | Output mains adapter 12Vdc/1000mA |

Security of memorized data

Unlimited, independent of battery charge conditions

Time

| | |
|---------------|----------------------|
| Date and time | Real time schedule |
| Accuracy | 1min/month max error |

Measured values storage - model HD2156.2

| | |
|------------------|---|
| Type | 2000 pages containing 10 samples each |
| Quantity | 20,000 sets of three measurements composed of pH or mV, χ , Ω or TDS or salinity and temperature. |
| Storage interval | 1s, 5s, 10s, 15s, 30s, 1min, 2min, 5min, 10min, 15min, 20min, 30min and 1h. |

Serial interface RS232C

| | |
|---------------------|--|
| Type | RS232C electrically isolated |
| Baud rate | Can be set from 1200 to 38400 bauds |
| Data bit | 8 |
| Parity | None |
| Stop bit | 1 |
| Flow Control | Xon/Xoff |
| Serial cable length | Max 15m |
| Print interval | Immediate or selectable between: 1s, 5s, 10s, 15s, 30s, 1min, 2min, 5min, 10min, 15min, 20min, 30min and 1h. |

USB interface - model HD2156.2

| | |
|------|---------------------------------|
| Type | 1.1 - 2.0 electrically isolated |
|------|---------------------------------|

Connections

| | |
|------------------------------------|---------------------------------------|
| pH/mV input | Female BNC connector |
| Conductivity and Temperature input | 8-pole male DIN45326 connector |
| Serial RS232C interface | 8-pole MiniDin connector |
| USB interface | MiniUSB B-type connector |
| Mains adapter | 2-pole connector (positive at centre) |

Measurement of pH by Instrument

| | |
|---|--|
| Measurement range | -2.000...+19.999pH |
| Resolution | 0.01 or 0.001pH selectable from menu |
| Accuracy | ± 0.001pH ±1 digit |
| Input impedance | > 10 ¹² Ω |
| Calibration error @25°C | 10offset>20mV Slope < 50mV/pH or Slope > 63mV/pH Sensitivity < 85% or Sensitivity > 106.5% |
| Temperature compensation automatic/manual | -50...+150°C |

Measurement of mV by Instrument

| | |
|--------------------|---------------------|
| Measurement range | -1999.9...+1999.9mV |
| Resolution | 0.1mV |
| Accuracy | ±0.1mV ±1digit |
| Drift after 1 year | 0.5mV/year |

| Measurement of conductivity | | Resolution | Measurement of total dissolved solids (with coefficient χ /TDS=0.5) | | Resolution |
|-----------------------------|--------------------------|------------------|--|-------------------------|------------|
| Measuring range Kcell=0.01 | 0.000...1.999 μ S/cm | 0.001 μ S/cm | Measuring range Kcell=0.01 | 0.000...19.999mg/l | 0.005mg/l |
| Measuring range Kcell=0.1 | 0.00...19.99 μ S/cm | 0.01 μ S/cm | Measuring range Kcell=0.1 | 0.00...19.99mg/l | 0.05mg/l |
| Measuring range Kcell=1 | 0.0...199.9 μ S/cm | 0.1 μ S/cm | Measuring range Kcell=1 | 0.0...199.9mg/l | 0.5mg/l |
| | 200...1999 μ S/cm | 1 μ S/cm | | 200...1999mg/l | 1mg/l |
| | 2.00...19.99mS/cm | 0.01mS/cm | | 2.00...19.99g/l | 0.01g/l |
| | 20.0...199.9mS/cm | 0.1mS/cm | | 20.0...99.9g/l | 0.1g/l |
| Measuring range Kcell=10 | 200...1999mS/cm | 1mS/cm | Measuring range Kcell=10 | 100...999g/l | 1g/l |
| Accuracy (conductivity) | $\pm 0.5\% \pm 1$ digit | | Accuracy (total dissolved solids) | $\pm 0.5\% \pm 1$ digit | |

| Measurement of instrument's resistivity | | Resolution | Measurement of salinity | | Resolution | |
|---|-----------------------------|--------------------------|-------------------------------------|--------------------------|--|--|
| Measuring range Kcell=0.01 | up to 1G Ω -cm (*) | | Measurement range | 0.000...1.999g/l | 1mg/l | |
| Measuring range Kcell=0.1 | up to 100M Ω -cm (*) | | | 2.00...19.99g/l | 10mg/l | |
| Measuring range Kcell=1 | 5.0...199.9 Ω -cm | 0.1 Ω -cm | Accuracy (salinity) | 20.0...199.9g/l | 0.1g/l | |
| | 200...999 Ω -cm | 1 Ω -cm | | $\pm 0.5\% \pm 1$ digit | | |
| | 1.00k...19.99k Ω -cm | 0.01k Ω -cm | | | | |
| | 20.0k...99.9k Ω -cm | 0.1k Ω -cm | | | | |
| | Measuring range Kcell=10 | 100k...999k Ω -cm | 1k Ω -cm | Temperature compensation | 0...100°C with α_T selectable from 0.00 to 4.00%/°C | |
| | | 1...10M Ω -cm | 1M Ω -cm | automatic/manual | | |
| Accuracy (resistivity) | $\pm 0.5\% \pm 1$ digit | | Reference temperature | 20°C or 25°C | | |
| | | | χ / TDS Conversion factor | 0.4...0.8 | | |
| | | | Cell constant K (cm ⁻¹) | 0.01, 0.7, 1.0 and 10.0 | | |

(*) The resistivity measurement is obtained from the reciprocal of the conductivity measurement: the indication of the resistivity, in the vicinity of the full scale, appears as in the following table

| K cell = 0.01 cm ⁻¹ | | K cell = 0.1 cm ⁻¹ | |
|--------------------------------|------------------------------|-------------------------------|------------------------------|
| Conductivity (μ S/cm) | Resistivity (M Ω -cm) | Conductivity (μ S/cm) | Resistivity (M Ω -cm) |
| 0.001 μ S/cm | 1000 M Ω -cm | 0.01 μ S/cm | 100 M Ω -cm |
| 0.002 μ S/cm | 500 M Ω -cm | 0.02 μ S/cm | 50 M Ω -cm |
| 0.003 μ S/cm | 333 M Ω -cm | 0.03 μ S/cm | 33 M Ω -cm |
| 0.004 μ S/cm | 250 M Ω -cm | 0.04 μ S/cm | 25 M Ω -cm |

Standard solutions automatically detected @25°C

- 147 μ S/cm
- 1413 μ S/cm
- 12880 μ S/cm
- 111800 μ S/cm

TECHNICAL DATA OF PROBES EQUIPPED WITH INSTRUMENT

2 and 4 electrode conductivity probes

| ORDER CODE | MEASUREMENT RANGE | DIMENSIONS |
|---|---|------------|
| SP06T | K=0.7 5 μ S...200mS/cm 0...90°C 4-electrode cell in Pohan/Platinum Max pressure 5bar | |
| SPT 401.001 not suitable for HD 2306.0 | K=0.01 0.04...20 μ S/cm 0...120°C 2-electrode cell AISI 316 - Teflon Max pressure 5bar | |
| SPT01G | K=0.1 0.1 μ S...500 μ S/cm 0...80°C 2-electrode cell in Glass/Platinum Max pressure 5bar | |
| SPT1G | K=1 10 μ S...10mS/cm 0...80°C 2-electrode cell in Glass/Platinum Max pressure 5bar | |
| SPT10G | K=10 500 μ S...200mS/cm 0...80°C 2-electrode cell in Glass/Platinum Max pressure 5bar | |

Temperature measurement of the instrument

| | |
|------------------------|-----------------|
| Pt100 measuring range | -50...+200°C |
| Pt1000 measuring range | -50...+200°C |
| Resolution | 0.1°C |
| Accuracy | ±0.1°C ±1 digit |
| Drift after 1 year | 0.1°C/year |

TECHNICAL DATA OF PROBES AND MODULES EQUIPPED WITH INSTRUMENT
Temperature probes Pt100 sensor with SICRAM module

| Model | Type | Application field | Accuracy |
|--------------------------------|-----------------------------|-------------------|---|
| TP472I | Immersion | -196°C...+500°C | ±0.25°C (-196°C...+300°C) ±0.5°C (+300°C...+500°C) |
| TP472I.0 1/3 DIN Thin Film | Immersion | -50°C...+300°C | ±0.25°C (-50°C...+300°C) |
| TP473PI | Penetration | -50°C...+400°C | ±0.25°C (-50°C...+300°C) ±0.5°C (+300°C...+400°C) |
| TP473P.0 1/3 DIN Thin Film | Penetration | -50°C...+300°C | ±0.25°C (-50°C...+300°C) |
| TP474C.I | Contact | -50°C...+400°C | ±0.3°C (-50°C...+300°C) ±0.5°C (+300°C...+400°C) |
| TP474C.0 1/3 DIN Thin Film | Contact | -50°C...+300°C | ±0.3°C (-50°C...+300°C) |
| TP475A.0 1/3 DIN Thin Film | Air | -50°C...+250°C | ±0.3°C (-50°C...+250°C) |
| TP472I.5 | Penetration | -50°C...+400°C | ±0.3°C (-50°C...+300°C) ±0.6°C (+300°C...+400°C) |
| TP472I.10 | Penetration | -50°C...+400°C | ±0.30°C (-50°C...+300°C) ±0.6°C (+300°C...+400°C) |
| TP49A.0 Class A Thin Film | Immersion | -70°C...+250°C | ±0.3°C (-70°C...-50°C) ±0.25°C (-50°C...+250°C) |
| TP49AC.0 Class A Thin Film | Contact | -70°C...+250°C | ±0.3°C (-70°C...-50°C) ±0.25°C (-50°C...+250°C) |
| TP49AP.0 Class A Thin Film | Penetration | -70°C...+250°C | ±0.3°C (-70°C...-50°C) ±0.25°C (-50°C...+250°C) |
| TP875.I | Globe-thermometer Ø150mm | -30°C...+120°C | ±0.25°C |
| TP876.I | Globe-thermometer Ø50mm | -30°C...+120°C | ±0.25°C |
| TP87.0 1/3 DIN Thin Film | Immersion | -50°C...+200°C | ±0.25°C |
| TP878.0 1/3 DIN Thin Film | Photovoltaic | +4°C...+85°C | ±0.25°C |
| TP878.1.0 1/3 DIN Thin Film | | | |
| TP879.0 1/3 DIN Thin Film | Compost | -20°C...+120°C | ±0.25°C |

Common characteristics

Temperature drift @ 20°C 0.003%/°C

4 wires Pt100 and 2 wires Pt1000 Probes

| Model | Type | Application field | Accuracy |
|----------------------------------|----------------|-------------------|----------|
| TP47.100.0 1/3 DIN Thin Film | 4 wires Pt100 | -50...+250°C | 1/3 DIN |
| TP47.1000.0 1/3 DIN Thin Film | 2 wires Pt1000 | -50...+250°C | 1/3 DIN |
| TP87.100.0 1/3 DIN Thin Film | 4 wires Pt100 | -50...+200°C | 1/3 DIN |
| TP87.1000.0 1/3 DIN Thin Film | 2 wires Pt1000 | -50...+200°C | 1/3 DIN |

Common features

Temperature drift @20°C

Pt100 0.003%/°C

Pt1000 0.005%/°C

ORDER CODES

HD2156.1: The kit is supplied with: instrument HD2156.1, 4 1.5V alkaline batteries, operating manual, case and DeltaLog9 software.

HD2156.2: The kit is supplied with: instrument HD2156.2 **datalogger**, 4 1.5V alkaline batteries, operating manual, case and DeltaLog9 software.

pH/mV probes, conductivity probes, temperature probes, standard calibration solutions for various types of measures, connection cables for pH electrodes with S7 connector, cables for data transfer to PC or printer have to be ordered separately.

HD2110CSNM: 8-pole connection cable MiniDin - Sub D 9-pole female for RS232C.

C.206: Serial connection cable with USB connector for PC and 8-pole MiniDin male connector for the instrument HD2156.1.

CP23: USB 2.0 connection cable type A - MiniUSB type B (not suitable for HD2156.1).

DeltaLog9: Software for download and management of the data on PC using Windows operating systems.

SWD10: Stabilized power supply 100-240 Vac/12Vdc-1A mains voltage


HD40.1: 24-column portable thermal printer, serial interface, 57mm paper width, four NiMH 1.2V rechargeable batteries, SWD10 power supply, instruction manual, 5 thermal paper rolls. It uses the cable HD2110CSNM (optional).

RCT: The kit includes 4 thermal paper rolls 57mm wide and 32mm in diameter.

BAT-40: Spare battery pack for HD40.1 printer with built-in temperature sensor.

HD22.2: Laboratory electrode holder composed of base plate with built-in magnetic stirrer, shaft and replaceable electrode holder. Suitable diameter 12mm. Powered by bench-top meters of the series HD22...with cable HD22.2.1 (optional) or power supplier SWD10 (optional).

HD22.3: Laboratory electrode holder composed of base plate. Flexible arm for free positioning. Suitable for electrodes with diameter 12mm.

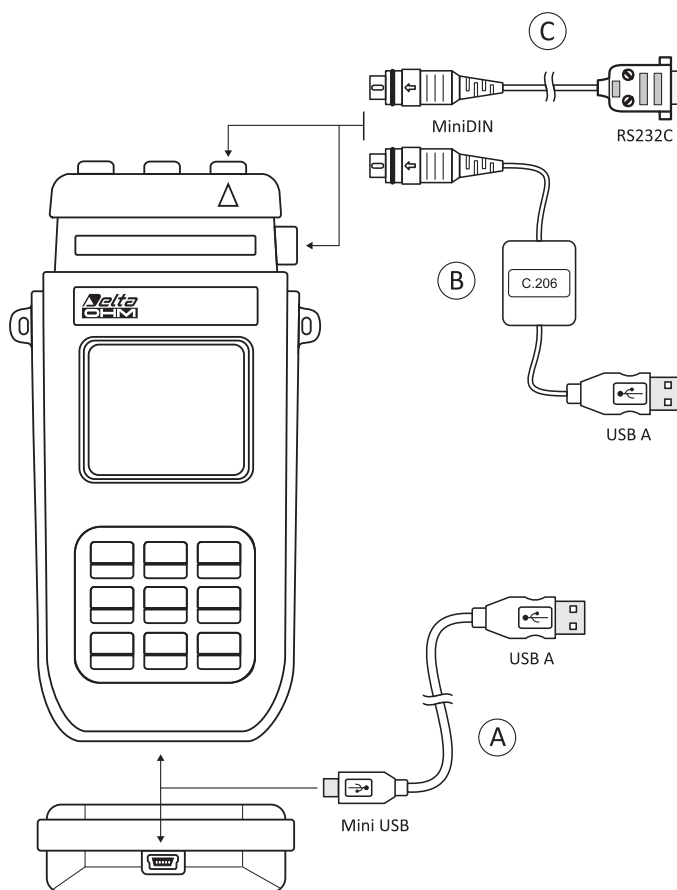

A For the models of portable data logger series **HD21XX.2** has been implemented with a new serial port miniUSB type HID (Human Interface Device).

When making the connection to the PC by the USB cable Type A - Mini USB B-type coded CP23, **no USB driver installation is requested.**

B For the connection of the models **HD21XX.1** to the RS232 port of your PC, the USB/serial converter is available (**code C.206**). The converter is equipped with its own drivers that have to be installed **before** connecting the converter to the PC (please see the details in the CD Rom supplied with the converter).

C The port with the MiniDIN connector which is present on every model is an RS232C type. By means of the cable coded HD2110CSNM, an RS232 port of a PC or the HD40.1. printer can be connected.

pH Electrodes

- KP 20:** Gel pH filled combined electrode for general use, with S7 screw connector, EPOXY body.
- KP 30:** Gel pH combined electrode for general use, 1m cable with BNC, EPOXY body .
- KP 50:** Gel pH combined electrode, porous Teflon ring junction, suitable for emulsions, demineralised water, waste water with S7 screw connector, glass body.
- KP 61:** 3 diaphragm liquid filled pH combined electrode for wine, milk, cream, etc., S7 screw connector, liquid reference filling, glass body.
- KP 62:** 1 diaphragm gel pH combined electrode for general use, pure water, varnishes, gel filled, S7 screw connector, glass body.
- KP 63:** liquid filled pH combined electrode for general use, varnishes, 1m cable with BNC, glass body.
- KP 64:** Liquid filled pH combined electrode, Teflon ring diaphragm, for wine, varnishes, emulsions, S7 screw connector, glass body.
- KP 70:** Pointed gel combined pH microelectrode diam. 6 x L=70 mm., with S7 screw connector, EPOXY body, glass tip, open junction for meat and cheese.
- KP 80:** Pointed gel pH combined electrode, with S7 screw connector, glass body, for cream, milk, viscous material, open junction.
- KP100:** Flat membrane gel combined pH electrode with S7 screw connector, glass body, for skin, leather, paper.

Characteristics and dimensions of the probes at page WA-76

- CP:** 1.5m extension cable with BNC/S7 connector for electrode without cable, thread S7.
- CP 5:** 5m extension cable with BNC/S7 connector for electrode without cable, thread S7.
- CP 10:** 10m extension cable with BNC/S7 connector for electrode without cable, thread S7.
- CP 15:** 15m extension cable with BNC/S7 connector for electrode without cable, thread S7.
- CE :** S7 screw connector for pH electrode.
- BNC:** female BNC for extension cable

ORP Electrodes

- KP 90:** REDOX PLATINUM liquid filled electrode with S7 screw connector, glass body.
- KP 91:** Gel REDOX PLATINUM electrode, 1m cable with BNC, EPOXY body for general purpose light duty.

Characteristics and dimensions of the probes at page WA-76

pH Buffer solutions

- HD8642:** Buffer solution 4.01pH - 200cc.
- HD8672:** Buffer solution 6.86pH - 200cc.
- HD8692:** Buffer solution 9.18pH - 200cc.

Redox Buffer solutions

- HDR220:** Redox buffer solution 220mV 0.5 l.
- HDR468:** Redox buffer solution 468mV 0.5 l.

Electrolyte solutions

- KCL3M** Ready to use solution for electrode refilling – 100 cc

Cleaning and maintenance

- HD62PT:** Diaphragm cleaning (tiourea in HCl) - 500ml.
- HD62PP:** Protein cleaning (pepsin in HCl) - 500ml.
- HD62RF:** Regeneration (fluorhydric acid) - 100ml.
- HD62SC:** Solution for electrode preservation - 200ml.

Conductivity probes

Please refer to the purchasing codes reported on the table of the probes on page WA-9.

Conductivity buffer solutions

- HD 8747:** Calibration solution 0.001 mol/l corresponding to 147 µS/cm at 25°C, 200cc.
- HD 8712:** Calibration solution 0.1 mol/l corresponding to 12,880 µS/cm at 25°C, 200cc.
- HD 8714:** Calibration solution 0.01 mol/l corresponding to 1413 µS/cm at 25°C, 200cc.
- HD 87111:** Calibration solution 1 mol/l corresponding to 111800 µS/cm at 25°C, 200cc.

Temperature probes equipped with SICRAM module

- TP472I:** Wire wound Pt100 sensor, immersion probe. Stem Ø 3 mm, length 300 mm. Cable length 2 m.
- TP472I.0:** Thin film Pt100 sensor, immersion probe. Stem Ø 3 mm, length 230 mm. Cable length 2 m.
- TP473P.I:** Wire wound Pt100 sensor, penetration probe. Stem Ø 4mm, length 150 mm. Cable length 2 m.

- TP473P.0:** Thin film Pt100 sensor, penetration probe. Stem Ø 4mm, length 150 mm. Cable length 2 m.
- TP474C.I:** Wire wound Pt100 sensor, contact probe. Stem Ø 4mm, length 230mm, contact surface Ø 5mm. Cable length 2 m.
- TP474C.0:** Thin film Pt100 sensor, contact probe. Stem Ø 4mm, length 230mm, contact surface Ø 5mm. Cable length 2 m.
- TP475A.0:** Thin film Pt100 sensor, air probe. Stem Ø 4mm, length 230mm. Cable length 2 m.
- TP472I.5:** Thin film Pt100 sensor, penetration probe. Stem Ø 6mm, length 500 mm. Cable length 2 m.
- TP472I.10:** Thin film Pt100 sensor, penetration probe. Stem Ø 6mm, length 1000mm. Cable length 2 m.
- TP49A.0:** Thin film Pt100 sensor, immersion probe. Stem Ø 2,7mm, length 150mm. Cable length 2 m. Aluminium handle
- TP49AC.0:** Thin film Pt100 sensor, contact probe. Stem Ø 4mm, length 150mm. Cable length 2 m. Aluminium handle
- TP49AP.0:** Thin film Pt100 sensor, penetration probe. Stem Ø 2,7mm, length 150mm. Cable length 2 m. Aluminium handle
- TP875.I:** Wire wound Pt100 sensor, 150mm diameter globe-thermometer equipped with handle. Cable length 2 m.
- TP876.I:** Wire wound Pt100 sensor, 50mm diameter globe-thermometer equipped with handle. Cable length 2 m.
- TP87.0:** Thin film Pt100 sensor, immersion probe. Stem Ø 3 mm, length 70 mm. Cable length 2 m.
- TP878.0:** Thin film Pt100 sensor, contact probe for solar panels. Cable length 2 m.
- TP878.1.0:** Thin film Pt100 sensor, contact probe for solar panels. Cable length 5 m.
- TP879.0:** Thin film Pt100 sensor, penetration probe for compost. Stem Ø 8 mm, length 1000 mm. Cable length 2 m.

Temperature probes without SICRAM module

- TP47.100.0:** Thin film Pt100 sensor, immersion probe. Stem Ø 3mm, length 230mm. Connection cable 4 wires with connector, length 2 m.
- TP47.1000.0:** Thin film Pt1000 sensor, immersion probe. Probe's Stem Ø 3mm, length 230mm. Connection cable 4 wires with connector, length 2 m.
- TP47:** Connector for Pt100 4-wire and Pt1000 2-wire probes without SICRAM module.
- TP87.100.0:** Thin film Pt100 sensor, immersion probe. Stem Ø 3mm, length 70mm. 4-wires connection cable with connector, length 1 m.
- TP87.1000.0:** Thin film Pt1000 sensor, immersion probe. Stem Ø 3mm, length 70mm. 2-wires connection cable with connector, length 1 m.

